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
Weegtechniek B.V.

uweshop.eu

KS

(Handleiding / Handbuch / Manual)

NL

D

GB

NL

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Thank you for choosing this RIBA scale.
designed for simple and straightforward use.

If you have any queries or experience any problems not addressed in the operating instructions, please contact your RIBA professional service partner.

Before take the scale into use, please read with care the information given in the operating instructions.
They contain important instructions for installation, proper use and maintenance of the appliance.

The manufacturer shall not be liable for damages arising out of failure to heed the following instructions:

When using electrical components under increased safety requirements, always comply with the appropriate regulations. Never perform work on the appliance while the power is switched on.

Inproper installation will render the warranty null and void.

The device is designed for use within buildings. Observe the permissible ambient temperatures for use (Technical information).

The device meets the requirements for electromagnetic compatibility. Do not exceed the maximum values specified in the applicable standards.

If you have any problems, contact your local RIBA professional service partner.

It is sufficient to clean the scale with a moist cloth and a regular cleaning agent. Do not use any scourers.

The scale does not require any routine maintenance. However, we recommend to check the scale's accuracy at regular intervals. The regularity of these checks is dependent on the level of use and the state of the scale. If any inaccuracies occur, please contact your local dealer or RIBA professional service partner.

Interference suppression of the unit corresponds to the valid directive.
The display value may be influenced by extreme electro-magnetic interference, e.g. operation of a radio unit or a mobile phone in the immediate vicinity of the unit. When the electrostatic interference is no longer present. The unit again be used for its intended purpose. You may need to switch it on again. If there is permanent electrostatic interference, we recommend earthing the platform.
The unit is a measuring instrument. Air currents, vibrations, rapid changes in temperature and direct sunlight may have an effect on the accuracy of the scale. The scale is protected to IP 54 protection class. Therefore, avoid high levels of humidity, vapours, aggressive liquid media and strong contamination.

Installation:

GB

Please check if you did receive following products:

* Balance

* Manual

Contact your supplier if you didn't receive some products.

Installation of the balance:

* Place the balance with the display faced towards you.

Specification:

Model:	KS
Weighing range:	6 kg
Division	1 gram
Platter size:	157x157 mm
Power:	Adapter / batteries
Operating temperature:	0°C - 40°C

Caution!

Please handle your balance with care.

Take notice of the following recommendations.

- * Liquids (like water) which come in contact with the balance can harm the accuracy.
 - * Overload of the balance can harm the accuracy of the balance.
 - * Extreme temperature changes can harm the accuracy of the balance.
 - * A clash with a heavy machine can make the balance useless.
 - * Changes in the balance may only be done by your supplier.
 - * We recommend you to let the balance check by your supplier every year.
 - * Only trained employees are allowed to use the balance.
 - * If you warm up the balance for 15 minutes before using, the accuracy of the balance will be better.
-
- * Always use the parts of your supplier.

Keyboard & Display:

- 1 UNIT KEY
- 2 ON/OFF KEY
- 3 ZERO KEY

Display unit:

Calibrate the scale

Only calibrate this scale if you are familiar with the technic and have the right aproved massweights

- 1 Switch off the scale
- 2 Press UNIT+ON/OFF keys together
Display shows the AD value
- 3 Press ON/OFF key
- 4 Put the displayed weight on the platter, confirm with On/Off key
- 5 PASS will appear and scale goes to weighingmode

Use tare function

Put the scale on a stable surface

Switch on the scale and wait for zero

put an empty container on the platter and press ZERO, display shows zero again

Your next weight will be NET weight

to erease tare weight just empty the platter and press ZERO again

Backlight and autom. Off settings

Press ZERO+ON/OFF together

Display shows b On, press ZERO to set backlite AUTO / or B OFF

Press UNIT keys to select P OFF, press ZERO to select P030 or P240 (power off setting)

Confirm the settings with UNIT key

Error messages

0_Id	Overload
Lo	Empty battery
Unst	Not stable
Err E	Memory error
Outz	Zero is out of range

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Maintenance:

Please disconnect the device from the operating voltage before cleaning.

Only use a cloth dampened with mild suds and not aggressive cleaning agents (solvents or similar).

Please ensure that fluids are not able to get into the device and rub off using a clean, soft cloth.

Loose sample residue/powder can be removed carefully using a brush or hand vacuum cleaner.

The operating company shall ensure that load-bearing components are checked by a competent person (your supplier) at maximum intervals of one year.